上海源叶生物科技有限公司
Shanghai yuanye Bio-Technology Co., Ltd
电话: 021-61312973 传真: 021-55068248
网址: www.shyuanye.com
邮箱: shyyssw@sina.com

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Description	ALC-0159 是一个聚乙二醇的脂质结合物, 可用作疫苗赋形剂。
In Vitro	Preparation of Lipid Nanoparticles Here we provide lipid molar ratios for LNPs in FDA-approved BNT162b2 (a COVID-19 mRNA vaccine). The molar ratio of lipids in this formulation is ALC-0315 : DSPC : Cholesterol : ALC-0159 = 46.3 : 9.4 : 42.7 : 1.6, and RNA to lipid weight ratio is 0.05 (wt/wt)[1]. A. Lipid Mixture Preparation 1. Dissolve lipids in ethanol and prepare 10 mg/mL stock solutions. The lipid stock solutions can be stored at 20°C for later use. Note 1: The ionizable lipid is usually a liquid. Due to the viscosity, it should always be weighed rather than relying on the autopipette volume. Note 2: Cholesterol in solution should be kept warm (>37°C) to maintain fluidity. Transfer the cholesterol solution promptly to avoid cooling. 2. Prepare the lipid mixture solution as described. For each mL of lipid mixture add the following: 560 µL of 10mg/mL ALC-0315, 261 µL of 10mg/mL Cholesterol, 117 µL of 10mg/mL DSPC, and 62 µL of 10mg/mL ALC-0159. Mix the solutions thoroughly to achieve a clear solution. This mixture contains 10 mg of total lipid. Note 3: The choice of lipids and ratios may be changed as desired and this will affect the LNP properties (size, polydispersity, and efficacy) and the amount of mRNA required. B. mRNA Preparation 1. Prepare a 166.7 µg/mL mRNA solution with 100 mM pH 5 sodium acetate buffer. Note 4: The lipid:mRNA weight ratio influences the encapsulation efficiency. Other weight ratios may be prepared as alternative formulations and should be adjusted accordingly by user. C. Mixing There are three commonly used methods to achieve rapid mixing of the solutions: the pipette mixing method, the vortex mixing method, and the microfluidic mixing method. All these mixing methods can be used for various applications. It is important to note that pipette mixing method and vortex mixing method may yield more heterogeneous LNPs with lower encapsulation efficiencies and is prone to variability. Microfluidic devices enable rapid mixing in a highly controllable, reproducible manner that achieves homogeneous LNPs and high encapsulation efficiency. Within these devices, the ethanolic lipid mixture and aqueous solution are rapidly combined in individual streams. LNPs are formed as the two streams mix and are then collected into a single collection tube. 1. Pipette Mixing Method: 1.1. Pipette 3 mL of the mRNA solution and quickly add it into 1 mL of the lipid mixture solution (A 1:3 ratio of ethanolic lipid mixture to aqueous buffer is generally used.) Pipette up and down rapidly for 20-30 seconds. 1.2. Incubate the resulting solution at room temperature for up to 15 minutes. 1.3. After mixing, the LNPs were dialyzed against PBS (pH 7.4) for 2 h, sterile filtered using 0.2 µm filters, and stored at 4°C. 2. Vortex Mixing Method:

上海源叶生物科技有限公司
Shanghai yuanye Bio-Technology Co., Ltd
电话: 021-61312973 传真: 021-55068248
网址: www.shyuanye.com
邮箱: shyyssw@sina.com

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